Centrifuge Operator's Manual
C-14000

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WARNING:

Use universal precautions when handling laboratory specimens. All human specimens of blood, blood compounds, or bodily fluids are potentially biohazardous and may contain Hepatitis B virus, Hepatitis C virus, Human immunodeficiency virus, or other disease-causing agents. For the safety of both the operator and service personnel, care should be taken when handling substances that are known to be toxic, radioactive, or contaminated with pathogenic microorganisms when using this centrifuge. When Risk Group II materials are used (as identified in the World Health Organization "Laboratory Bio-Safety Manual") use of a Bio-Seal should be employed. More than one level of protection must be provided in the case of materials of a higher group. The use of flammable or explosive materials, as well as those materials which chemically react vigorously, is prohibited. Any use of the equipment in a manner not specified in these instructions may impair the level of protection provided by the equipment.

Model Description

The Centrifuge is a continuous-duty, electronically controlled, variable-speed, centrifuge equipped with a lid safety interlock system. It offers the user control over many parameters, including variable runtime, speed control, and deceleration. The exclusive rotor allows for quick and easy sample loading and complete horizontal separation. Samples can be viewed safely through the transparent lid. Entry into the centrifuge is restricted during operation by the safety interlock system. Electronic push buttons and a backlit display easily control centrifugation parameters.

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Intended Use

This centrifuge is an IVD accessory, and therefore subject to the former EU IVD Directive 98/79/EC and the current EU IVD Regulation 2017/746. It is a laboratory product designed to separate components by the generation of relative centrifugal force. It separates human and animal samples – such as blood, urine, and other body fluids - collected in an appropriate specimen receptacle, either alone or with reagents or other additives. As a general-purpose laboratory centrifuge, it is designed to also run other containers filled with chemicals (non-flammable, non-explosive, non-volatile, and non-highly reactive only), environmental samples, and other non-human body samples. This centrifuge should be operated by trained personnel only. Any use other than those intended by the Manufacturer is explicitly prohibited. The maximum sample density is 1.15 grams/mL.

Supplied Equipment*

- One (1) 8-Carrier Fixed Angle Rotor
- Eight (8) 22 mL Inserts
- Eight (8) 11 mL Inserts
- Eight (8) Caps
- One (1) Operator's Manual
- One (1) Power Cord

*The rotor and rotor accessories are rated for a rotational frequency of 4,000 RPM.

The use of any cord other than what is supplied by the manufacturer may not carry an adequate rating and is therefore prohibited.

Features

- Incorporates a unique tube holder, insert, and cap
- A lid safety switch that prevents the centrifuge from operating unless the lid is closed and latched
- Locking lid that allows entry into the centrifuge only after the rotor has completely stopped
- Clear lid for safe observation of samples
- Electronically controlled operation
- Push-button operation

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Overall Dimensions (H x W x D):

370mm x 330mm x 175mm

Timer: Electronic, 15 seconds to 99 minutes, preset to 10 minutes

Voltage Requirements: Frequency: 115/ 230 VAC (+/- 10%)

Permitted Environmental Conditions

Ambient Temperature During Operation:

60°F - 90°F (16°C - 32°C)

Maximum Relative Air Humidity: 90%

Setup Location

Unpack the centrifuge and verify that all of the supplied equipment is present. Choose a setup location that meets the following criteria:

- A benchtop clearance height of 20" is required to open the lid.
- The clearance envelope is the space around the centrifuge that is required for safety.
- Choose a setup location that will allow for a clearance envelope of at least 24" x 24", (with the centrifuge at the center).
- Do not position the centrifuge in such a way that it would be difficult to disconnect the power in case of emergency.
- No person or hazardous material shall be permitted in the clearance envelope during operation.
- The operator time within the envelope shall be limited to the time necessary for loading, unloading, and centrifuge operation only.
- Proper ventilation is necessary to prevent the overheating of samples as well as premature failure of the centrifuge.
- Choose an area that will allow unencumbered airflow.
- No adjustment is necessary for leveling the centrifuge; however, the surface should be flat and level.
- Be sure the outlet is always within reach as the line cord is the means of emergency disconnection!

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Initial Setup Procedure

If any problems are found during the initial setup procedure, refer to the troubleshooting section.

Plug the female end of the supplied line cord into the power input module located on the rear of the centrifuge. Plug the male end into an approved electrical outlet. For electrical safety, the unit must always be properly grounded.

Flip the switch on the power input module located on the rear of the centrifuge to the ON (I) position. For safety purposes, the locking system is always activated. To deactivate the system, (to insert or retrieve samples), press the OPEN button on the control panel.

Spin the rotor by hand; check for free and level rotation. If the rotor does not spin freely, refer to the troubleshooting section.

Start a centrifugation cycle by pushing the START button. The unit will accelerate to full speed.

Listen to the sound of the centrifuge. A smooth whirring sound should be heard. If there are any loud or unusual sounds, stop the centrifuge by pushing the STOP button immediately and refer to the troubleshooting section.

Push the STOP button. The motor should slow to a stop.

- The lid should remain locked until the rotor has nearly stopped. Once the rotor has stopped, the interlock system will become disengaged.

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Operation

NOTE: Follow the initial setup procedure.
Push the OPEN button and then open the lid.

If using 11/12ml tubes, place tubes in the white inserts and place the inserts in the black tube holders. If using 22/23ml tubes, place in the black tube holders and load the tube holders into the rotor. Place a cap over each tube holder containing a sample tube or balance tube. Be sure to follow the rules for balanced loads below.

Close the lid. If the lid is not completely latched, the centrifuge will not operate!

1. Turn on the machine by pushing the START button on the control panel.
2. The centrifuge should begin to spin.
3. After time has elapsed, the rotor will slow to a complete stop.
4. The locking mechanism will disengage allowing entry into the rotor chamber. If it does not, refer to the troubleshooting section.
- 5.
6. Remove the samples.

Balanced Loads

Your centrifuge must contain a balanced load to work properly. Use the following rules when loading the rotor. Spinning balanced loads will extend the life of the machine and produce better results.

Opposing tube holders must be identical
Opposing tube holders must be empty or loaded with equally weighted samples.

If an odd number of samples is to be spun, fill a balance tube with water to match the weight of the unpaired sample and place it across from this sample.

Place a cap on the black tube holder containing the balance tube and the sample tube to balance the load.

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Care and Preventative Maintenance

With proper care and maintenance, your centrifuge will provide years of laboratory service. For proper care, the following steps should be taken:

Provide Adequate Ventilation: For cooling purposes, the centrifuge draws in ambient air through the air intake cover on the top of the lid and exhausts this air in the rear of the base. The centrifuge should be placed on a hard smooth surface for good air circulation.

Always Spin Balanced Loads: Make certain that you are always spinning a balanced load. The centrifuge has a unique counter-balanced motor mounting design which, along with its rubber suction feet, produces excellent vibration dampening. However, out-of-balance loads may break glass test tubes and may produce unsatisfactory separation results. Proper load balancing will improve sample separation and extend the life of the centrifuge. Refer to the section on balanced loads for additional information on balancing the load.

Keep the Tube Holders Clean: NOTE: Always follow the safety guidelines of your laboratory to properly clean up and/or dispose of materials if a substance known to be potentially toxic, radioactive, or contaminated with a pathogenic microorganism is spilling in or on the centrifuge. Small glass fragments left in the tube holder after a tube breakage may adhere to the next test tube inserted in that holder. When this tube is handled, these fragments may puncture protective gloves and lacerate the operator's fingers or hands. The remaining fragments may provide stress points on subsequent tubes and result in additional breakage. If a tube breakage occurs, carefully remove the tube holder. Properly dispose of the sample and tube fragments and thoroughly clean both the inside and outside of the tube holder. Replace the tube holder in the rotor.

Motor and Electrical Maintenance: The centrifuge uses a brushless permanent split capacitor AC motor. It should not need servicing for the life of the centrifuge. The electrical components are selected for high reliability and should not need service.

Remove Accessories Before Moving: All tube holders, samples, and caps must be removed from

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the rotor chamber before transporting or storing the centrifuge to prevent damage and injury.

Cleaning and Disinfection

To prolong the life of the centrifuge cleaning and disinfection is recommended every six months, or whenever there is a spillage or tube breakage. Contaminants must be removed immediately, or corrosion and premature degradation of components can occur.

Unplug the centrifuge before cleaning.

Apply cleaning solutions with a towel or cloth. Do not submerge the centrifuge in water or other cleaning solutions as this will cause damage and void the warranty.

ONLY isopropyl alcohol, soap, and water, or a 10% (5500 PPM) bleach solution should be used for cleaning and disinfection of the centrifuge and accessories.

All surfaces must be dried immediately after cleaning and disinfecting.

TBQ Germicidal products should not be used, as they will cause damage to the centrifuge and void the warranty.

The use of fully/partially halogenated hydrocarbons, ketones, esters, ethers, benzyls, ethyl benzenes, and all other chemicals not prescribed by the manufacturer shall not be used as they may cause damage to the rotor chamber, rotor, tube holders, accessories, and centrifuge exterior and void the warranty.

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Troubleshooting

Problem: The rotor does not spin freely.

Solutions:

- Make sure nothing has fallen into the rotor chamber.
- If nothing is obstructing the rotor, contact Estar for further assistance.

Problem: Excessive noise when the machine is running.

Solutions:

- Check to see that the load is balanced.
- Make sure that nothing has fallen into the rotor chamber.
- Make sure that the nut in the center of the rotor is tight.

Problem: The centrifuge does not run.

Solutions:

- Check the electrical outlet.
- Make sure the lid latch is closed.

- Contact Estar for further assistance.

Transportation

During transportation, accessories can become dislodged from the rotor and fall into the rotor chamber where they can remain unnoticed. If the centrifuge needs to be transported to another site, make certain that there are no tube holders or accessories in the rotor or rotor chamber.

Repaired Units

For verification of the safe state of the centrifuge following factory service or repair, refer to the date on the Factory Calibration located on the back of the centrifuge. This is the date that the repaired centrifuge was the last factory tested and calibrated. For additional servicing information and technical support, contact Estar.

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Safety

Lid Safety Interlock System:

In addition to the Lid Safety Switch, The Centrifuge has a true "0 RPM" lid-locking system. The lid safety interlock system keeps the lid locked at all times, (even during a power failure), and requires that the rotor is at rest to unlock the lid. The centrifuge will not allow entry into the rotor chamber unless the centrifuge has power and the rotor is stopped. To open the lid, make sure that the centrifuge is plugged in and, with the rotor stopped, press the OPEN button.

Emergency Rotor Chamber Entry:

In the event of power failure, it may be impossible to unlock the lid by conventional means. In this case, entry into the rotor chamber may be made by disengaging the locking mechanism (located on the bottom under the front right of the machine). Push the mechanism to the right and open the lid. If the unit is damaged, contact Estar.

15, Hamerkava St., Holon
5885111, Israel
Tel: +972-3-5596414
Fax: +972-3-5596424
www.estar-medical.com

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